



## A NEW WAY to SAFELY DISPOSE OF Ethidium Bromide Solutions

A new product safely traps used solutions of ethidium bromide for disposal. Gravity fed, it can treat 3-5 liters of EtBr waste ( $0.5 \mu\text{g/ml}$ ) to a minimum detectable limit of  $0.025 \mu\text{g/ml}$ .

Simply attached our cartridge, via the provided tubing, to the source of the waste and start it flowing by pulling on a syringe from the opposite end.

The cartridge will develop a reddish color as the EtBr begins to adsorb onto the bed and the color front will advance until all the EtBr has been absorbed. At that point, allow several milliliters of additional fluid to flow through the cartridge, disconnect it and properly dispose of it with other solid waste.

The EtBr is trapped on the cartridge and contact with lab personnel is minimized or eliminated.

ETBR	Ethidium Bromide Trap, Pk/1	\$30
ETBR-5	Ethidium Bromide Trap, Pk/5	120

*These prices have not been finalized and should be at or below those shown*